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光阻法 (Light Obscuration Particle Count Test)

For preparations supplied in containers with a nominal volume of more than 100 mL, apply the criteria of Test 1.A. For preparations supplied in containers with a nominal volume of less than 100 mL, apply the criteria of Test 1.B. For preparations supplied in containers with a nominal volume of 100 mL, apply the criteria of Test 1.B. [NOTE—Test 1.A is used in the Japanese Pharmacopoeia.] If the average number of particles exceeds the limits, test the preparation by the Microscopic Particle Count Test. Test 1.A (Solutions for parenteral infusion or solutions for injection supplied in containers with a nominal content of more than 100 mL)—The preparation complies with the test if the average number of particles present in the units tested does not exceed 25 per mL equal to or greater than 10 μm and does not exceed 3 per mL equal to or greater than 25 μm . Test 1.B (Solutions for parenteral infusion or solutions for injection supplied in containers with a nominal content of less than 100 mL). The preparation complies with the test if the average number of particles present in the units tested does not exceed 6000 per container equal to or greater than 10 μm and does not exceed 600 per container equal to or greater than 25 μm .

显微镜法 (Microscopic Particle Count Test)

For preparations supplied in containers with a nominal volume of more than 100 mL, apply the criteria of test 2.A. For preparations supplied in containers with a nominal volume of less than 100 mL, apply the criteria of test 2.B. For preparations supplied in containers with a nominal volume of 100 mL, apply the criteria of test 2.B. Test 2.A – Solutions for infusion or solutions for injection supplied in containers with a nominal content of more than 100 mL. The preparation complies with the test if the average number of particles present in the units tested does not exceed 12 per milliliter equal to or greater than 10 μm and does not exceed 2 per milliliter equal to or greater than 25 μm . Test 2.B – Solutions for infusion or solutions for injection supplied in containers with a nominal content of less than 100 mL. The preparation complies with the test if the average number of particles present in the units tested does not exceed 3000 per container equal to or greater than 10 μm and does not exceed 300 per container equal to or greater than 25 μm .

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